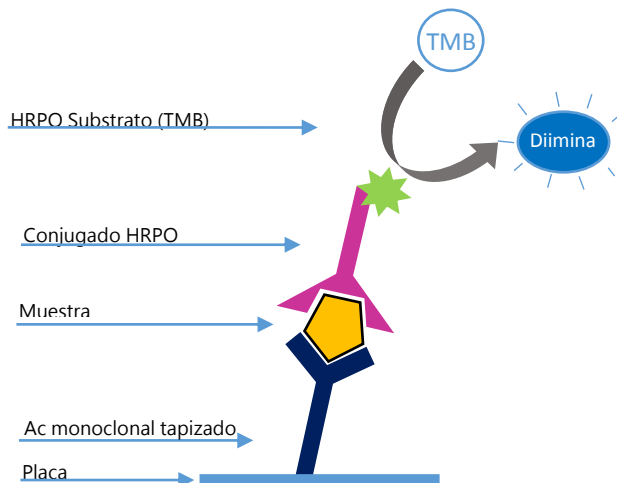
## INGEZIM CASEINA

R.30.CAS.K2

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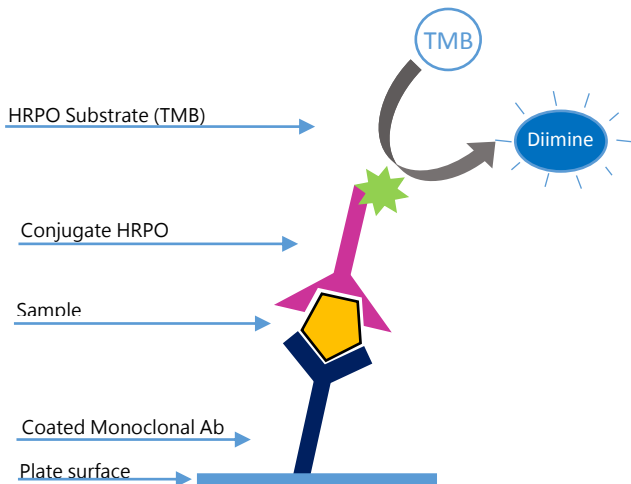
## INGEZIM CASEÍNA

R.30.CAS.K2

**Ingezim CASEINA** is an immunoenzymatic assay which uses specific monoclonal antibodies of casein from milk. **Limit of Detection: 0.25 ppm.**

### TECHNICAL BASIS OF THE KIT

1. Plates are supplied coated with Monoclonal Antibodies specific for CASEIN. Samples are added to the wells and incubated.
2. If the sample contains the antigen, it will bind to the antibody specific for casein coating the plates.
3. When MAb-PO specific of casein is added, it will bind to the antigen of the sample.
4. Binding is detected by a colorimetric reaction after the addition of the substrate

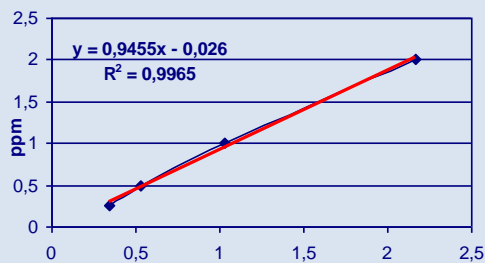
### APLICACION

This assay is a rapid and reliable test that significantly reduces the time required to screen food products and environmental samples for the presence of CASEIN residues.

Results can be read visually (qualitative interpretation), or using an ELISA reader (quantitative interpretation).

### SENSITIVITY

The assay is able to detect **0.25ppm** of CASEIN, by using a linear regression curve. It leads to more precise results (mainly at low concentrations).

Regression curve

### SPECIFICITY

The assay is specific for casein from cow, sheep and goat milk.

### VARIABILITY

Both the standard curve and an extracted sample were analyzed by duplicate, using 3 different lots. The average results, deviation and VC are showed in the table below.

Standards	Average	$\sigma$	vc
2 ppm	2,05	0,03	1,42
1 ppm	1,1	0,02	1,81
0,5 ppm	0,63	0,04	6,51
0,25 ppm	0,35	0,03	8,36
Sample	0,49	0,01	2,24

### KIT COMPOSITION

- Microtitration plate of 48 wells
- Vial with Negative Control
- CASEIN Standards (4)
- Vial with PO-Conjugate
- Bottle with Washing Solution
- Bottle with substrate (TMB)
- Bottle with stop solution
- Bottle with Extraction Solution

PRODUCT MANUFACTURED BY INGENASA

IT-73840 IT-73780  
ISO 14001:2015 9191.INGE  
ISO 9001:2015 9175.ING2

SELF LIFE: **12 months**  
Stored at 4°C

02/02/2018