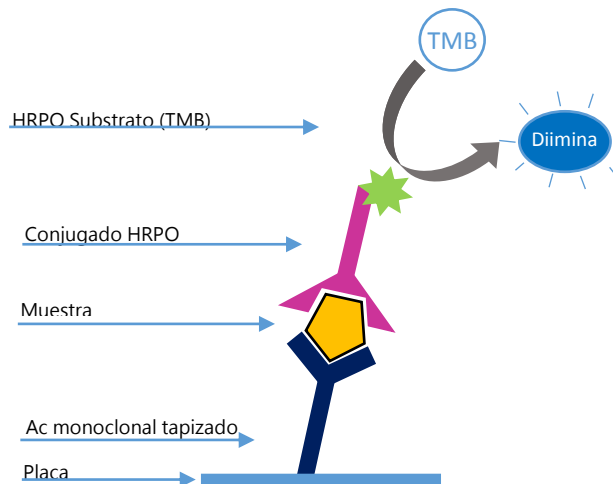
INGEZIM B-LACTOGLOBULINA

R.30.BLG.K2

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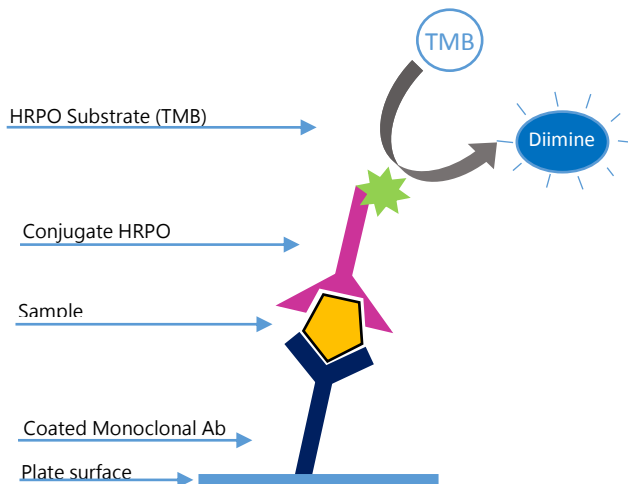
INGEZIM B-LACTOGLOBULINA

R.30.BLG.K2

Ingezim b-lactoglobulina is an immunoenzymatic assay which uses specific antibodies of b-lactoglobulin (BLG), the main protein in lacto serum. **Limit of Detection: 0.1 ppm.**

TECHNICAL BASIS OF THE KIT

1. Plates are supplied coated with Antibodies specific for BLG. Samples are added to the wells and incubated.
2. If the sample contains the antigen, it will bind to the antibody specific for BLG coating the plates.
3. When Ab-PO specific of BLG is added, it will bind to the antigen of the sample.
4. Binding is detected by a colorimetric reaction after the addition of the substrate

APLICACION

This assay is a rapid and reliable test that significantly reduces the time required to screen food products and environmental samples for the presence of BLG residues.

Results can be read visually (qualitative interpretation), or using an ELISA reader (quantitative interpretation).

SENSITIVITY

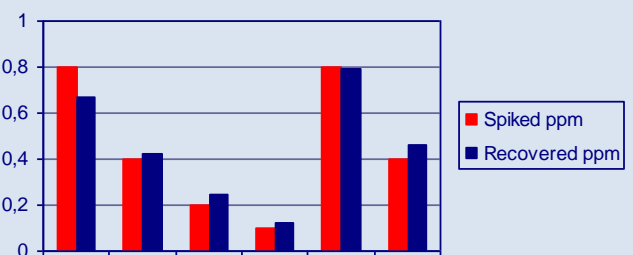
The assay is able to detect **0.1ppm** of b-lactoglobulin, by using a linear regression curve. It leads to more precise results (mainly at low concentrations).

Regression curve

RECOVERY DATA

Spike samples are prepared at different contamination level, and extracted several times. The average recovery data are showed below.

KIT COMPOSITION

- Microtitration plate of 48 wells
- Vial with Negative Control
- BLG Standards (4)
- Vial with PO-Conjugate
- Bottle with Washing Solution
- Bottle with substrate (TMB)
- Bottle with stop solution
- Bottle with Extraction Solution

PRODUCT MANUFACTURED BY INGENASA

UNDER PATENT LICENCE FROM CSIC AND HERO ESPAÑA

SELF LIFE: **12 months**
Stored at 4°C

02/02/2018