INgezim CTV R.20.CTV.K2

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

Ed.020217

Ingezim CTV R.20.CTV.K2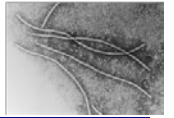

Ingezim CTV is a Double Antibody Sandwich immunoenzymatic assay for the detection of Citrus Tristeza Virus.

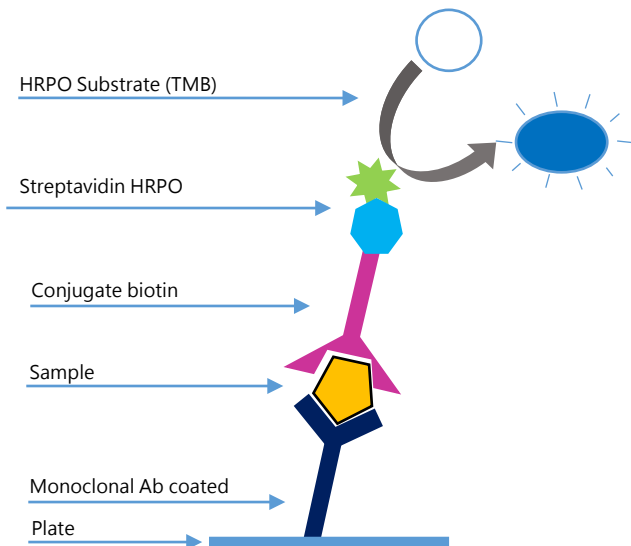TECHNICAL BASIS OF THE KIT

1. Plates are coated with monoclonal antibodies specific to CTV. Samples are added and incubated.
2. If the sample contains CTV it will bind to the specific MAb of CTV coating the plate.
3. After the addition of a biotin-conjugated MAb specific of CTV, it will bind to the antigen in case the sample was positive.
4. In order to detect this reaction, an Alkaline-Phosphatase conjugated Streptavidin is added.
5. The complex Streptavidin-biotin is detected by the development as a colorimetric reaction after the addition of an Alkaline-Phosphatase specific substrate.

APPLICATION

Detection of Citrus Tristeza Virus in citric plant extracts.

SPECIFICITY AND SENSITIVITY

1. Extracts of Mexican lime plants infected with 20 CTV isolates, which had previously been catalogued by 2 DAS ELISAs which use polyclonal antibodies (PAb-ELISA-DAS), were analyzed by Ingezim CTV. The controls used consisted of a mixture of young shoots from healthy Mexican lime (negative) and a mixture of young shoots from 350 Mexican lime seedlings infected with THE strain T-308. Ingezim CTV detected as positive all infected samples and the positive control. The healthy plants controls were negative, matching with the expected results.
2. Extracts from 1795 field samples were analyzed by Ingezim CTV in parallel with 2 PAb-ELISA_DAS. Ingezim CTV detected 584 samples as positive and 1199 as negative, showing 100% sensitivity and 99% specificity in relation to Pac-ELISA-DAS used.

Production and characterization of monoclonal antibodies specific for Citrus Tristeza Virus and their use for diagnosis. J. Gen. Virol. 1986;67, 91-96. C. Vela et al.

COMPOSITION OF THE KIT

- CTV specific monoclonal antibody for coating
- CTV specific Biotin conjugated monoclonal antibody
- Alkaline-Phosphatase- Streptavidine conjugate

1000 DETERMINATIONS **5000 DETERMINATIONS**

PRODUCT MANUFACTURED IN INGENASA

SHELF LIFE: 24 months
Stored at -20°C

Ed.020217